
Set content

1. Microscope

2. Flexible tripod arm with handle

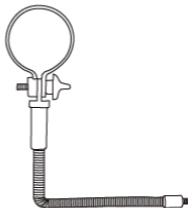3. Tripod base

4. MicroUSB cable


5. 4x anti-slipping pads


Device components


1. Lighting adjustment
2. Power/photo
3. Zoom and focus adjustment
4. Spacer sleeve with dust cover
5. Digital image magnification
6. White LED Backlight
7. Lens
8. Battery Charging Led
9. Power LED
10. Charging/Data microUSB socket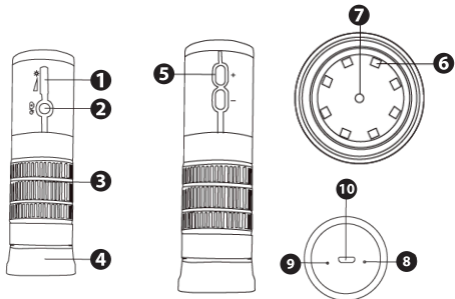


Assembling

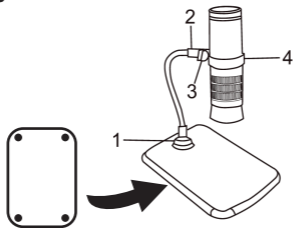
1.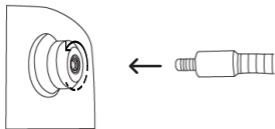


2.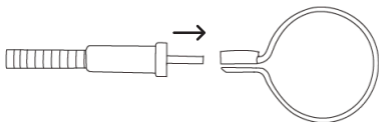

3.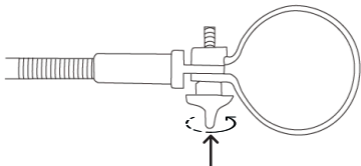

4.

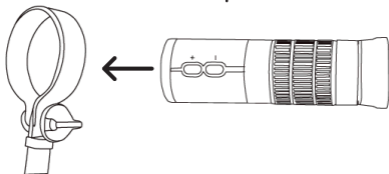
Software

The DM WIFI app is used to support MT4105 on IOS and Android mobile devices. Download it from store system and install:

How to use

1. Enable mt4105 with power button.
2. Search your smartphone for a WIFI network called DM WIFI and connect to it. You don't need a password.
3. If you receive the message „This network does not have internet access. Still want to connect to it?“, check the box „Don't ask for this network again“ and confirm „YES“
4. Open the DM WIFI app to get started.

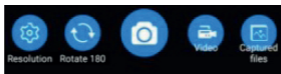5. Use the knob (3) to adjust the sharpness of the observed detail.

Note. It is possible to use the microscope in cooperation with your PC and the AMCAP shareware application available on the product page www.media-tech.eu

Battery charging

Charge with the included microUSB cable connected to the USB port on your computer or a standard USB 5V 1A charger

Note. Do not use fast QC3.0 chargers

When charging the battery, the charging light (9) is lighting. When the battery is fully charged, which lasts approx. hours, the LED will turn off.

Due to continuous development, the specifications and appearance of the device are subject to change without prior notice. The current version of the user manual and technical support available on www.media-tech.eu.

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GB Note on environmental protection:

After the implementation of the European Directive 2012/19/UE in the national legal system, the following applies: Electrical and electronic devices may not be disposed of with domestic waste. Consumers are obliged by law to return electrical and electronic devices at the end of their service lives to the public collecting points set up for this purpose or point of sale. Details to this are defined by the national law of the respective country. This symbol on the product, the instruction manual or the package indicates that a product is subject to these regulations. By recycling, reusing the materials or other forms of utilising old devices, you are making an important contribution to protecting our environment.

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