MICROSCOPE USB 500X

MT4096

- USB PC microscope (USB camera)
- Adjustable magnification 50-500x
- CMOS sensor
- Photos up to 30 Mpix (6324x4742ppi) after interpolation
- 300Kpix optical resolution
- Video recording: up to 25k / s at 640x480ppi
- Backlight: 8 white LEDs
- Light adjustment with a potentiometer
- High-quality metal tripod
- USB 2.0 interface
- Compatibility: Win 8.1 /10 32-64bit

User Manual

Thank you for purchasing our MT4096 MICROSCOPE USB 500X, it is a slim type tool and can be easily used in different fields including:

- Textile industrial for textile inspection
- Printing inspection
- Industrial inspection: PCB, precision machinery...
- Educational purpose
- Hair examination
- Skin examination
- Microbiological observation
- Jewelry & coin (Collections) inspection
- Visual Assistance

Notes before use

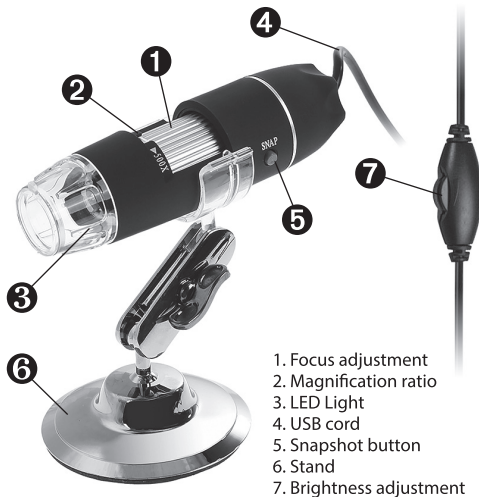
- Don't disassemble the digital Microscope or change the interior parts, it can cause damage.
- Don't clean the Microscope with alcohol organic solvents
- Don't touch the lens with your fingers.
- Avoid outdoor use if possible.
- Storage temperature, 0°C~40°C, Humidity: 45%RH~85RH%.
- In case the product gets wet, leave PC connection immediately. And do not disassemble or dry by hair dryer. Send to repair center if the digit microscope was effected by liquid or other elements.
- Measurement Data only for reference.

Product Outlook & Standard parts

Set completion

- Microscope with cable (1pcs)
- Stand (1pcs)
- CD ROM (Driver, Measurement software, User Manual)

Parts of microscope


1. Focus adjustment
2. Magnification ratio
3. LED Light
4. USB cord
5. Snapshot button
6. Stand
7. Brightness adjustment

How to install the Driver

U500X Digital Microscope can be installed after inserting attached CD (requires operation system above win2000 version), please make sure to plug in the device before installation.

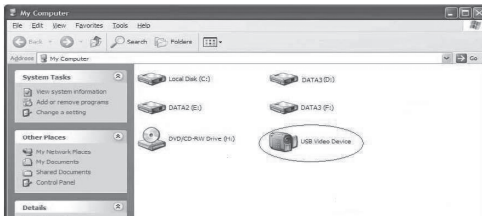1. Connect the Portable digital microscope by USB2.0 cable to your computer, double click 'my computer' icon on your computer desktop, below interface will be shown up.

2. Double click the **Video device icon** as above red marked in your computer. Now it is ready to use.

If you can't find the Video device, or if your computer can't recognize it, please follow below steps to install the driver. Insert the attached CD, system will auto run to the interface as below and follow the steps to complete the installation.

NOTE: If the computer can't auto run the disc, you need to operate it yourself.

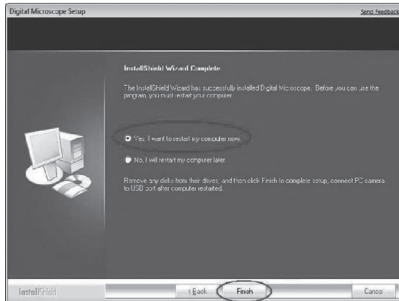
Driver installation

Click **Driver** to continue, the screen will show the installation progress.

Click "Next", as followings:

Select 'Yes, I want to restart my computer now', and click "finish".

After restart your computer, Plug-in the USB port of Digital Microscope into Computer USB port. Following the user guide by steps and complete by click "Next" -> "Continue anyway" -> "Finish".

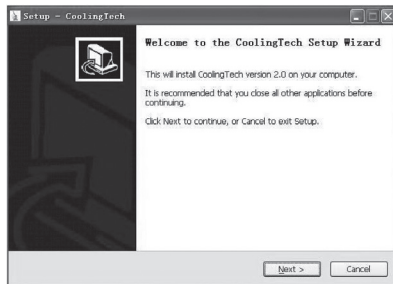
NOTE. Coolingtech software (Measurement) support Operation System Win7, Win8 but doesn't support Windows10. Please open "Driver" file folder in the CD and start the "setup.exe", you can use usb microscope in Windows10 after software "amcap" installed.

In Windows10, if happen error on operating "set camera size" or "capture picture" when using coolingtech measurement software, please do as the followings:


1. After installing the software, find "CoolingTech_Microscope" by the installation file folder path ,if you choose the default path, please select the file folder "CoolingTech_Microscope" under the file folder "C:\Program Files (x86)\CoolingTech";
2. Right click on the folder "CoolingTech_Microscope" and select "properties" -> "safe" -> "Edit" button, and then click"add (D)" button, please enter "everyone" under the blank of "enter the object name to choose.;" and then click"OK" button;
3. Please click "Everyone" under"group or user name (G):" , ->"Permission of Everyone" ->"complete control" ->"allow";
4. Click on the "application" and "OK" button ,then the coolingtech measurement software can be used normally.





Measurement installations

Click **Measurement** to continue, as followings:

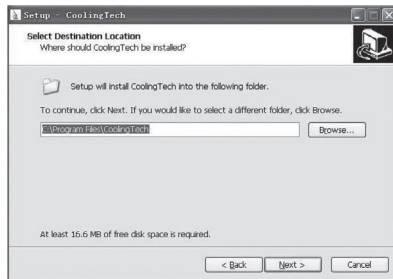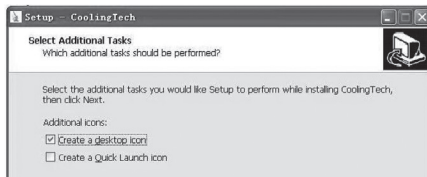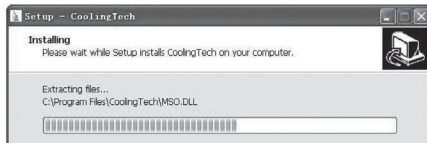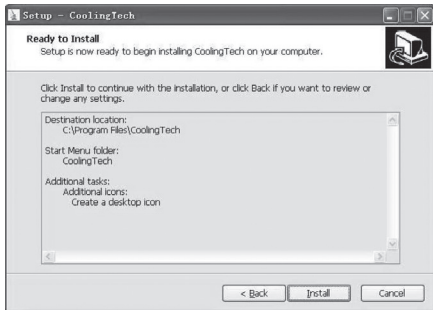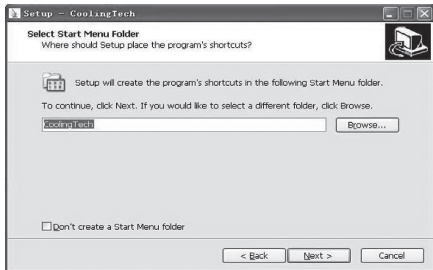Click "next" to continue, as followings.


Click "Finish" and the installation completes.


After the installation completes, double click icon at computer desktop to start using the software.

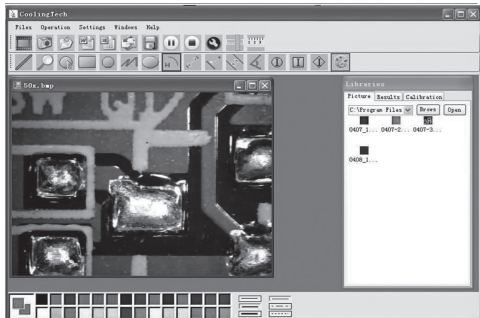Plug-in the devices, Display in operation window.


Click to open the operation system.

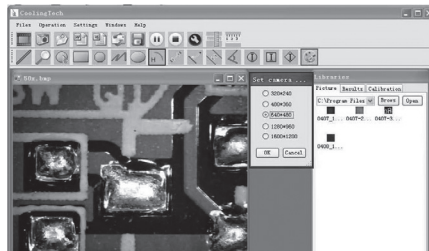Choose the Image resolution from 320*240 to 1600*1200, default at VGA Mode 640*480

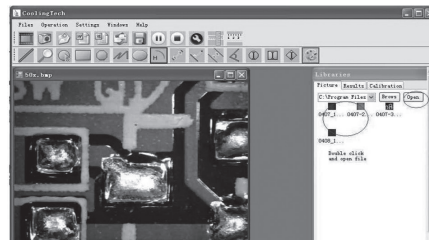Click to use measurement function.


Click to open pictures, and click to save pictures.

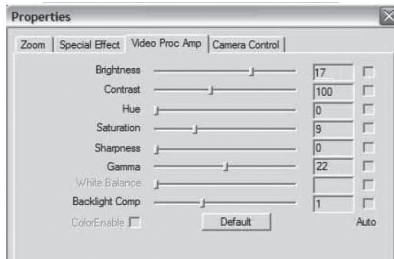Image Parameter including Image setting and special effects.

Video PLAY (Pause), STOP

Magnifier

Click Ruler

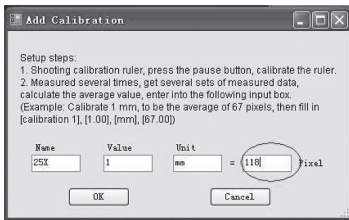
Click to, start calibration make sure certain accuracy.

Our experience told us that measurement could be controlled at 99% high precision. Measurement could be caused from human eye distortion and defocus. More practice is helpful to measurements.

To make the measurements of the captured images, you must adjust the calibration value yourself, using the attached calibration card.


Key in the Measure data (Pixel), Unit and Click'OK'


Warning: Please do calibration again once Magnification Ratio changed.

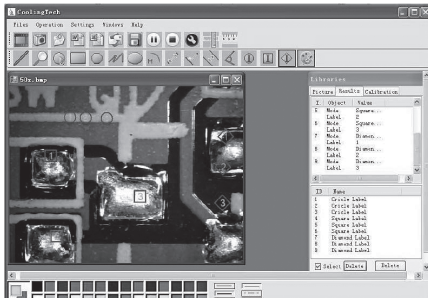Color setting of display character.


Click to export word or excel format.

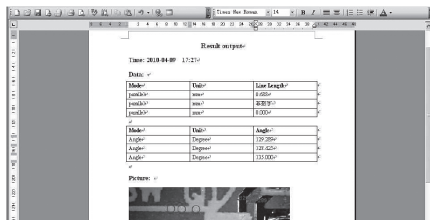
Measure: line, ring, square, polygon, set-line length, oval, arc, point to point distance, point to line distance, parallel line, angle.

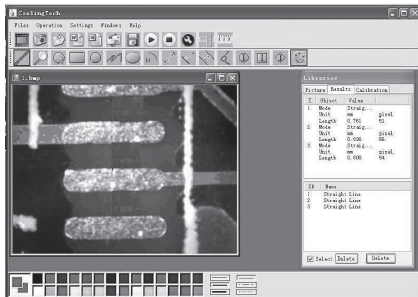
Line measurement:


Ring measurements:

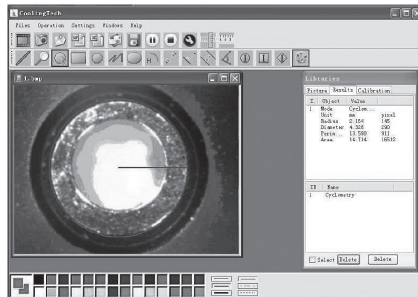Square measurement:

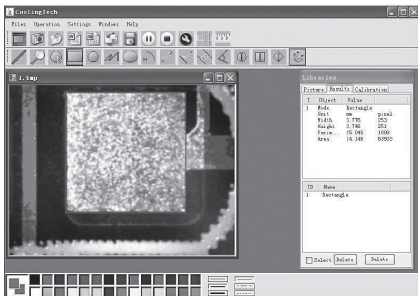Arc measurements:

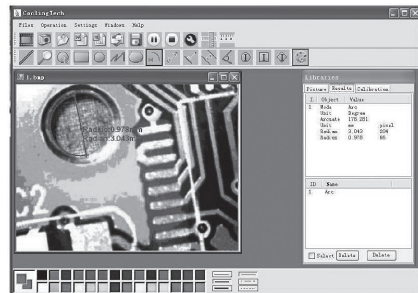*Due to continuous development specification and appearance of product are subject to change without prior notice.
For technical support please visit www.media-tech.eu.*

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EN Environmental guidance:

The device is marked with the symbol of a crossed out waste container, in accordance with the European Directive 2012/19/EC on waste electrical and electronic equipment. Products marked with this symbol should not be disposed of or disposed of after the end of their useful life together with other household waste. The user is obliged to dispose of waste electrical and electronic equipment by delivering it to a designated point where such hazardous waste is recycled. The collection of this type of waste in separate places and the proper process of their recovery contribute to the protection of natural resources. Proper recycling of waste electrical and electronic equipment has a beneficial effect on human health and the environment. In order to obtain information on where and how to dispose of waste electrical and electronic equipment in an environmentally safe way, the user should contact the relevant local authority, the waste collection point or the point of sale where he purchased the equipment.

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